


# Protein Demonstration Lab Sheet


1. Circle the sample(s) you predict contain protein.

apple juice    egg white    cooking oil    water    milk    lemon-lime soda

2. Add 5 mL of each sample to their own individual test tube. Record which sample is in each test tube and use colored pencils to show what the samples look like in each test tube.

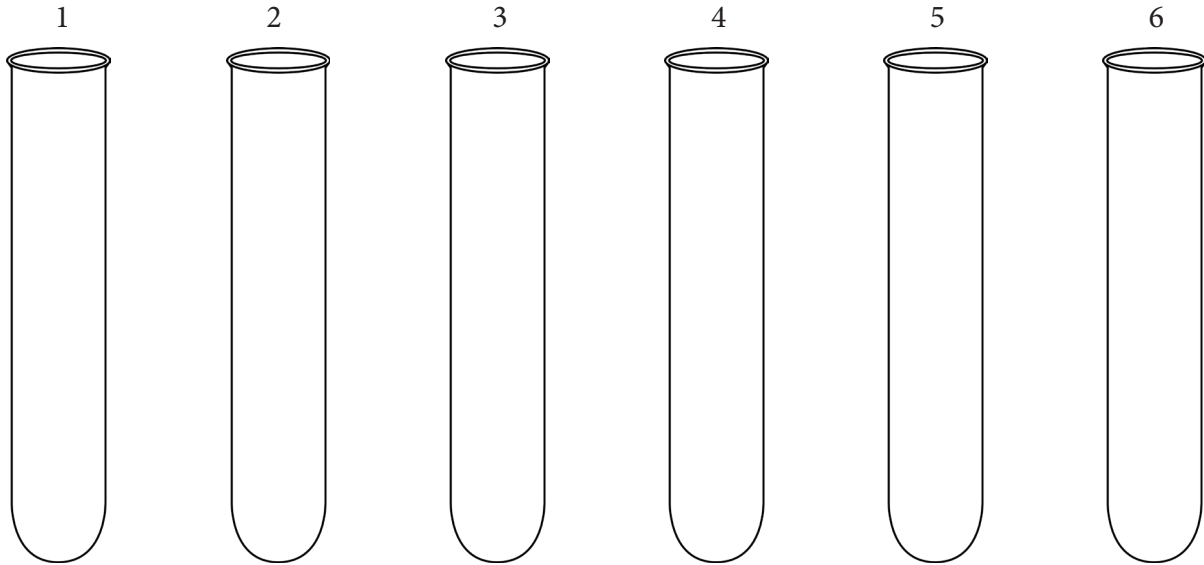

3. Add 15 drops of Biuret reagent to each test tube, stir well, and allow to sit for 5 minutes. Record your observations of what the solution in each test tube looks like after the Biuret reagent is added.

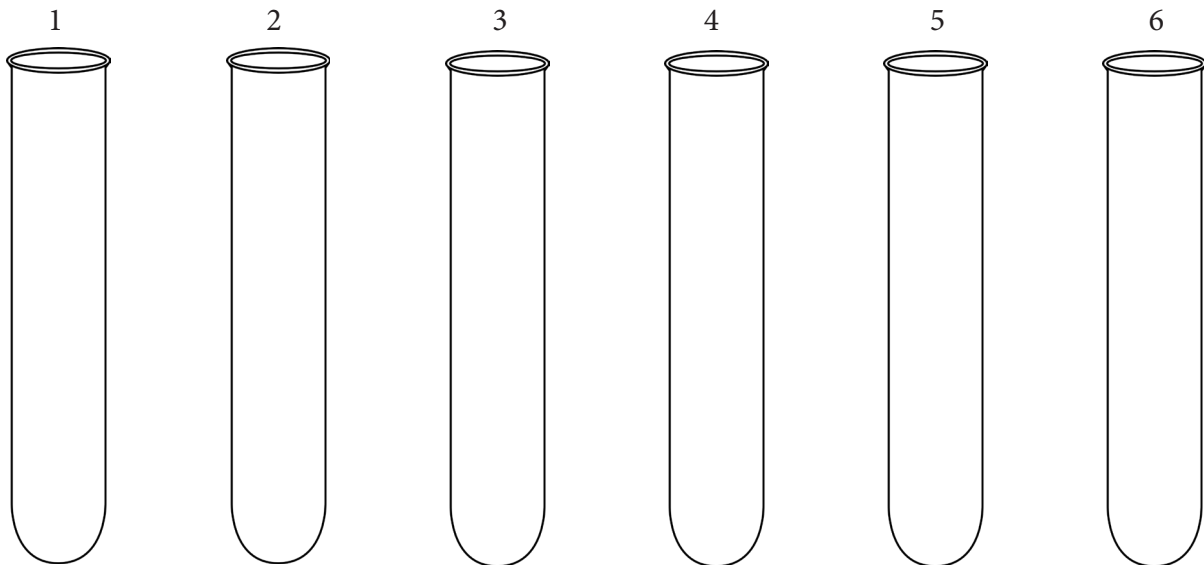

4. Which sample(s) contain protein? \_\_\_\_\_